# Installation Procedure of IRsolution

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**Read this manual thoroughly before you use the instrument.  
Keep this manual with care so that you can use it any time you need.**

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ANALYTICAL & MEASURING  
INSTRUMENTS DIVISION

Copyright © Shimadzu Corporation 2004. All rights are reserved, including those to reproduce this publication or parts there of in any form without permission in writing from Shimadzu corporation.

- Information in this publication is subject to change without notice and does not represent a commitment on the part of the vendor.
- Any errors or omissions which may have occurred in this publication will be corrected as soon as possible.
- Shimadzu does not have any responsibility for errors or injures resulting from following the instructions in this manual.

**Windows® is registered trademark of Microsoft Corporation.**

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**This is an instruction manual to install the IRsolution software which controls a FTIR and manipulate FTIR data. Refer to the Instruction manual – Operation guide of the IRPrestige-21/FTIR-8400S or Help message of the IRsolution software for software operation.**

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# Installation Procedure of IRsolution (English)

## Contents

### I Installation Procedure in English

- 1 Inspecting Standard Accessories .....2
- 2 Installation of Instrument Drivers.....3
  - 2.1 Installation of an IRPrestige-21 Driver program .....3
  - 2.2 Installation of an FTIR-8000 series Driver program .....3
- 3 Installation of IRsolution.....4
- 4 Installation of Supplemental Disk .....10
- 5 Selecting an Instrument .....11
- 6 User Management.....12
- 7 Uninstallation of the IRsolution .....13

# 1 Inspecting Standard Accessories

After unpacking, confirm that all parts are included in accordance with the standard accessory list shown below.

Table 1.1 Standard Accessories

	Part name	Quantity	Part No.
1	IRsolution software	1	206-72926-91
2	Instruction Manual IRPrestage-21/FTIR-8400 (Operation Guide (Japanese))	1	206-94966
3	Instruction Manual IRPrestage-21/FTIR-8400 (Operation Guide (English))	1	206-94967
4	Registration Card (Only for Japan)	1	206-72406-91
5	Guide to CLASS customer Service (Only for Japan)	1	223-03860
6	Certificate of Compliance	1	206-95059
7	Supplemental Disk	1	206-72924-91
8	Installation Procedure of IRsolution (This manual)	1	206-95080

## 2 Installation of Instrument Drivers

When your computer controls the IRPrestige-21/FTIR-8000 series, an instrument driver program must be installed. When the PC is for the data manipulation and/or printing, not for FTIR control, installation of instrument driver is not needed.

### 2.1 Installation of an IRPrestige-21 Driver program

A driver program for the IRPrestige-21 should be installed to control it.

1. Wire the IRPrestige-21 and your PC by refereeing to the section [Wiring the Cabled] of the Instruction Manual – User’s System Guide.
2. Turned them ON, and then Windows automatically recognizes the IRPrestige-21 with Plug & Play function.
3. Then Windows requests installation of the driver program. When Windows asks the place to find the driver, check [CD-ROM drive] in the Driver setup wizard dialog box.
4. Insert the [IRsolution] installation CD-ROM into your CD-ROM drive. When the Installation program of IRsolution starts, cancel it by clicking the <Cancel> button.
5. Follow the message on the screen. Then Windows automatically installs necessary drivers.
6. Install IRsolution software in next step.

#### Note

“FTIR-2010”, which is displayed on the found hardware wizard dialog box, is a module name of IRPrestige-21.

### 2.2 Installation of an FTIR-8000 series Driver program

A driver program for the FTIR-8000 series should be installed to control it.

1. Wire the FTIR-8000 series and your PC by refereeing to the section [Wiring the Cabled] of the Instruction Manual (User’s System Guide).
2. Turned the FTIR ON and then turn the PC ON. Windows automatically recognizes the FTIR-8000 series with Plug & Play function.
3. Then Windows requests installation of the driver program. When Windows asks the place to find the driver, check [CD-ROM drive] in the Driver setup wizard dialog box.
4. Insert the [IRsolution] installation CD-ROM into your CD-ROM drive. When the Installation program of IRsolution starts, cancel it by clicking the <Cancel> button.
5. Follow the message on the screen. Then Windows automatically installs necessary drivers.
6. Install IRsolution software in next step.

## 3 Installation of IRsolution

Install the IRsolution software with the following procedure.

1. When the IRsolution Install CD-ROM is in the CD-ROM drive, Click the <Start> button, and then select the [Run] command.

Select “setup.exe” on the CD-ROM drive, and then click <Run> button.

The Install program starts.

2. In another case, prepare the [IRsolution] install CD-ROM.

Turn on the power of the personal computer to start up the Windows.

If any software is running, terminate all of them.

Insert the CD-ROM into the CD-ROM drive.

The setup program automatically starts up. Execute “setup.exe” if setup program does not run automatically.

### Note

- If an old version of IRsolution is already installed, a message shown in Fig. 3.2 is displayed. Un-install old version of IRsolution first, and then install latest IRsolution again.

Fig. 3.1 A message when old version of IRsolution is installed

- English version of IRsolution install program is activated on another version of Windows than Japanese.
3. “Preparing Setup” dialog box and then “Welcome to the InstallShield Wizard for IRsolution” dialog box are displayed. Click the [Next] button.

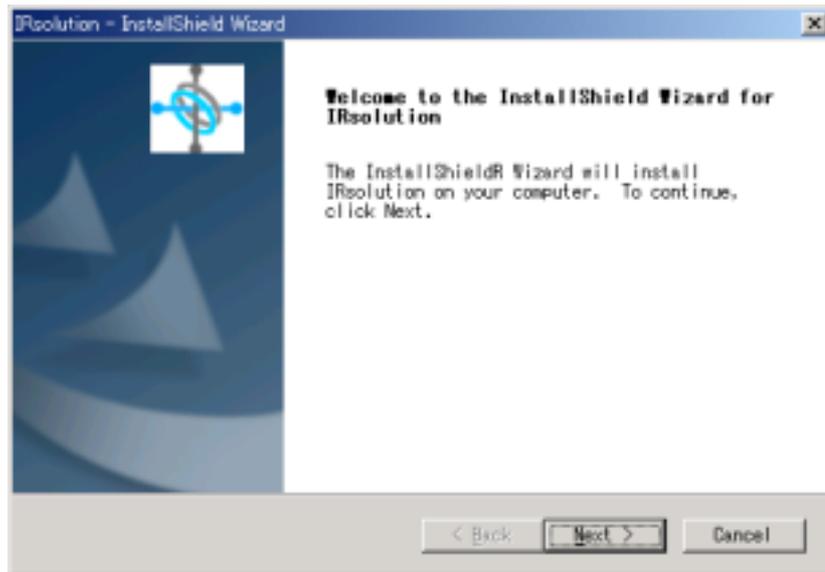
Fig. 3.2 “Welcome” dialog box

4. “Ready to Install the Program” dialog box is displayed. Click the [Install] button.

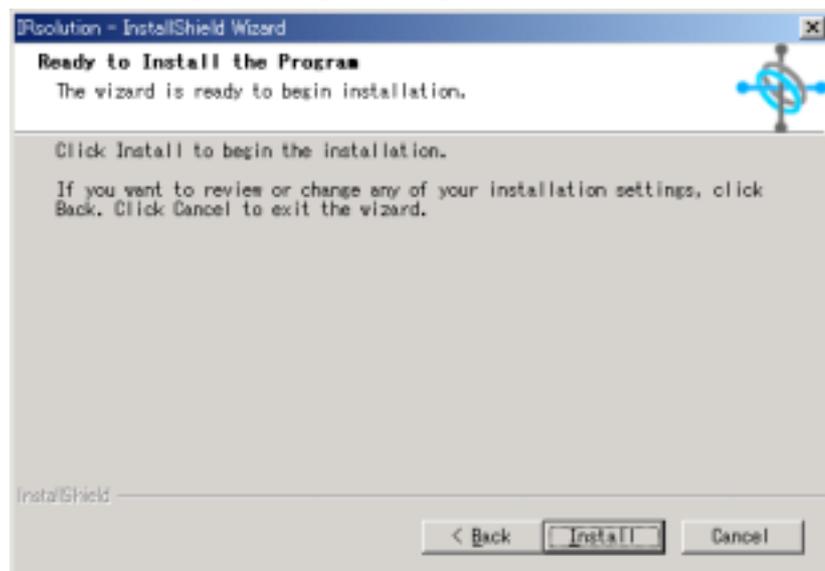
Fig. 3.3 “Ready to Install the Program” dialog box

5. Installation starts. Installation procedure is displayed on the “Setup Status” dialog box.
6. The [Please give in CD key for this product] window opens.

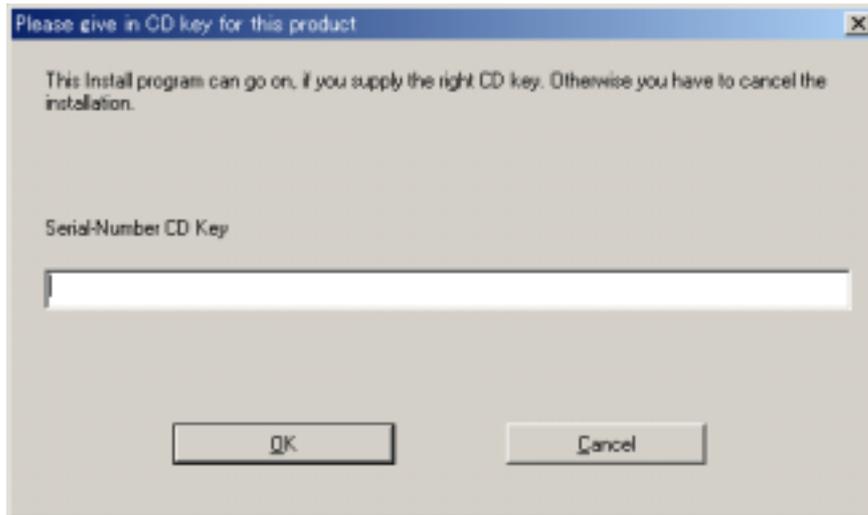
Fig. 3.4 Please give in CD key for this product

7. Input the CD Key written on the CD-ROM package onto the [Serial Number CD Key]. If input CD Key is not correct, you cannot proceed to next step. After inputting information, click the <OK> button.
8. Proceed following prompts on the screen. Then the [Chose Destination Location] window opens. Specify the destination folder to be installed. The default folder is used usually. Click the <Browse> button to change the destination.

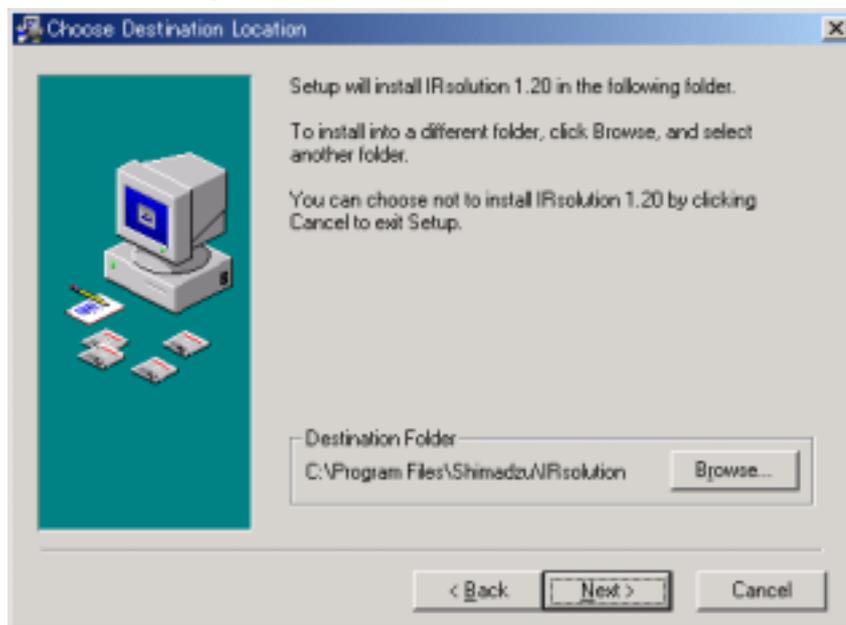
Fig. 3.5 Chose Destination Location

9. Click the <Next> button to proceed to next step, then the [Backup Replaced File] window opens. Click the <Next> button with the default setting.

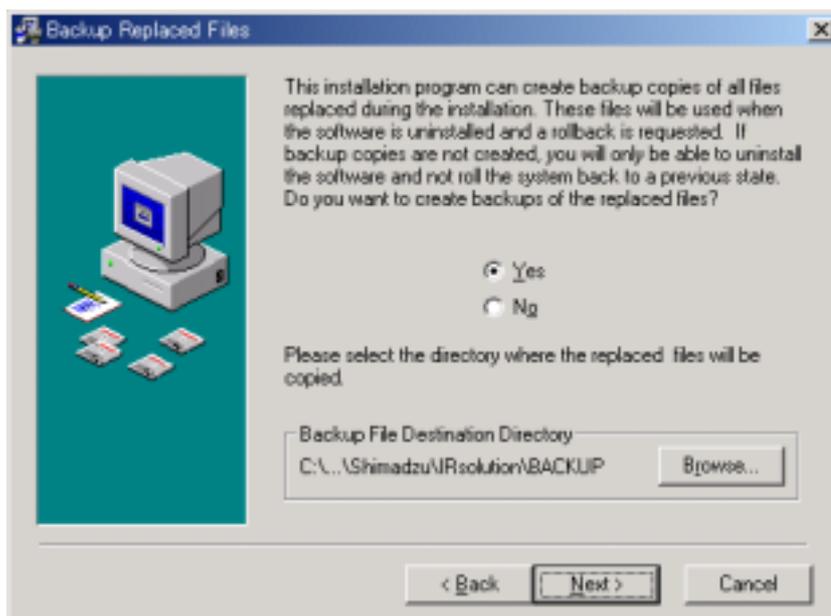
Fig. 3.6 Backup Replaced File

- The [Select Component] window opens Select components of the IRsolution main components to be installed. Following components can be installed.

Table 3.2 Setup Components (IRsolution)

Setup	Description
Program	The IRsolution main body
Help	Help message file of the IRsolution
GLP mode	Installs the IRsolution on the GLP mode to support the GLP/GMP. Electric Signature function is available on the GLP mode. You cannot switch the GLP/non-GLP mode after installation.
Support for hardlock protected libraries	Installs the hardlock driver to use protected libraries with a protection key (dongle).
DataExtractor support	IS checked, when used with the CLASS-Agent/IRsolution Agent.
Support for Sadtler hardlock Rainbow	Installs the hardlock driver to use Sadtler libraries

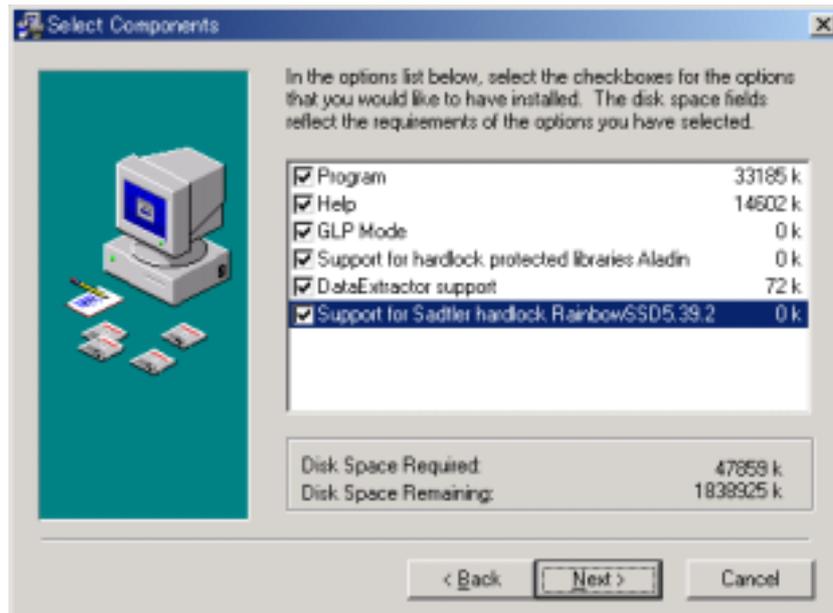
Fig. 3.7 Select Component

11. Put check marks on the necessary components. Click the <Next> button after setting.
12. The [Special Shortcuts] window opens. Put a check mark on the place you want to have shortcut icons. The default setting is usually used. Click the <Next> button.

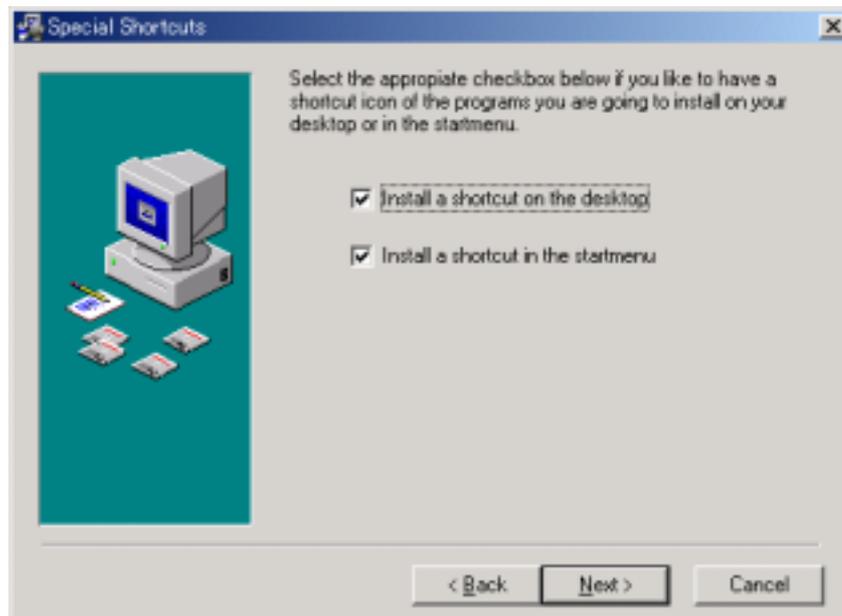
Fig. 3.8 Special Shortcuts

13. Proceed with following the prompts on the screen. The [Installation Completed] window opens when the IRsolution main body is installed correctly.
14. Click the <Finish> button to install the Alternation Check Program and Sample Files.
15. When all selected components are installed, the [Install Shield Wizard Complete] window opens. Click <Finish> button.

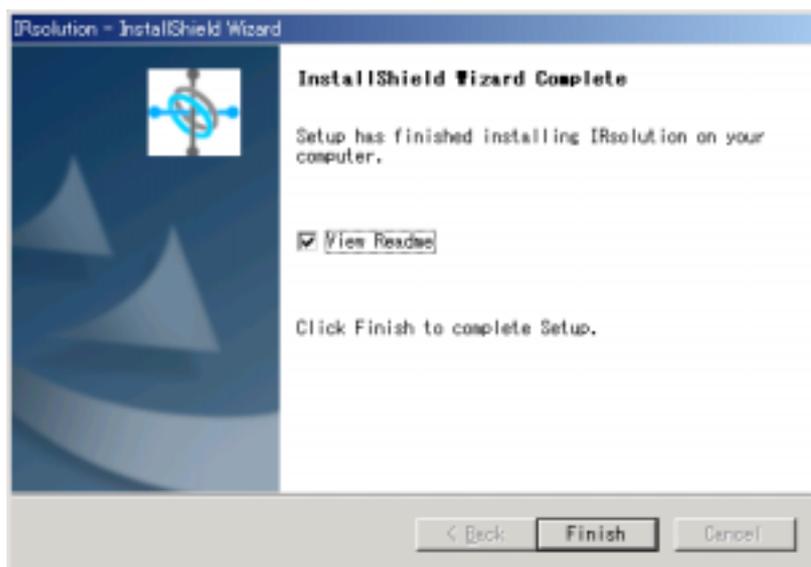
Fig. 3.9 Install Shield Wizard Complete

16. When a check mark is on the “View Readme”, Readme file is opened.
17. Install the Supplemental Disk in the next step.

Latest information for IRsolution software is described in ReadmeE.htm in both of Installation CD-ROM and Supplemental Disk. Please refer both of them before using IRsolution software.

**Caution**

This caution is related to the users who selects Windows XP as an operational system, and log on the IRsolution with privileges of Power Users group. In the described condition above, security settings of the folder are a little bit complicated and data files created by other personnel may not be usable. To avoid this problem, the installation program of IRsolution creates a data folder C:\Program Files\Shimadzu\IRsolution\data (where C represents the drive Windows was installed.) to which security settings are modified so that everyone can share data files. As it may not be possible to create folders or files, if they are created at the location other than C:\Program Files\Shimadzu\IRsolution\data and its sub folders, it is necessary to modify folder securities using tools provided by Windows.

## **4 Installation of Supplemental Disk**

Install the Supplemental Disk to update files and programs for the IRsolution software. Some version of IRsolution does not need installation of Supplemental Disk.

1. Insert the supplemental disk into disk drive in your computer.
2. Click the <Start> button, and then select [Run] command from the menu.
3. Input "(drive name):¥Setup.exe" in the file name and click <Run> button.
4. Run a set-up program, and updated files and programs are installed.

Latest information for IRsolution software is described in ReadmeE.htm in both of Installation CD-ROM and Supplemental Disk. Please refer both of them before using IRsolution software.

## 5 Selecting an Instrument

Setup the model of Instrument to be connected.

1. Activate the IRsolution software. Double click the IRsolution icon on the desktop or select the IRsolution menu on the Start menu.
2. Set up the model of instrument to connect with the IRsolution. Select the [Environment]-[Instrument Preferences]-[Instruments] menu before initializing the FTIR. Select [IRPrestige-21] as the Scanning FTIR Module. If you are using FTIR-8000 series, It's not necessary to change the setting.

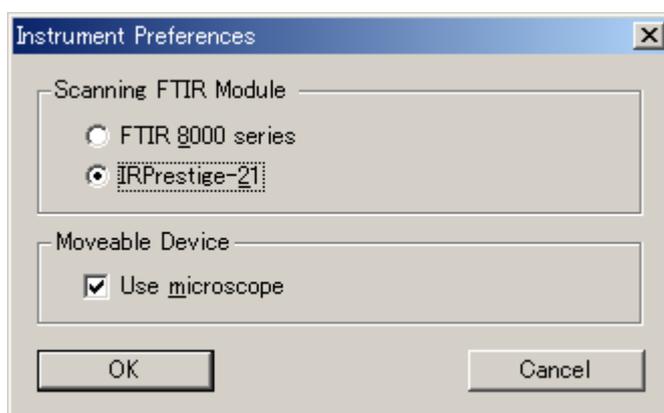
Fig. 5.2 [Instrument Preferences] Dialog box

FTIR 8000series	Connects to the FTIR-8000 series.
IRPrestige-21	Connects to the IRPrestige-21.

## 6 User Management

IRsolution has a security function by management of users with User ID and Password. Administrator should activate the security function by User ID and Password after installation of IRsolution by use of the [Admin]-[Security] command. Put a check mark on the “Password Restrictions” on the [Security] dialog box.

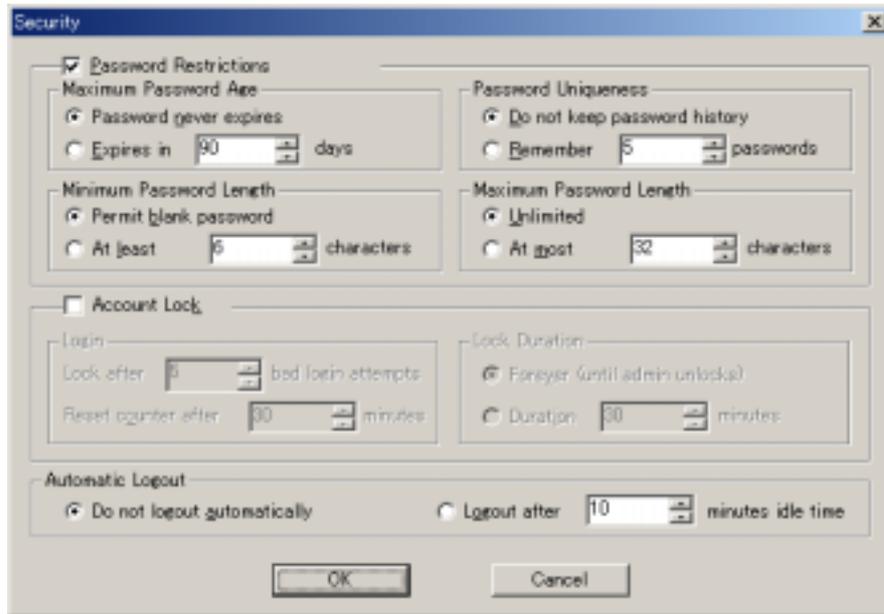
Fig. 6.1 [Security] dialog box

Clicking the <OK> button to open the message [There is no active user “xxxxxx”. The application has to be closed. Do you want to continue and close the application?]. Click the <Yes> button to close IRsolution once. The dialog box shown in Fig.6.2 is displayed when IRsolution is run again. Select “Administrator” as User name and nothing is input on the “Password” section, then click the <OK> button.

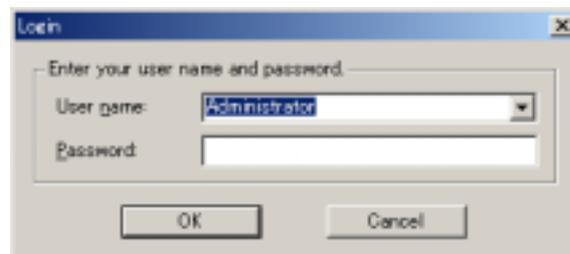
Fig. 6.2 [Log On] dialog box

Change the password for the Administrator by the [Admin]-[Security] command. Then manage the User and his/her rights of operations.

## 7 Uninstallation of the IRsolution

Follow the steps to uninstall the IRsolution software.

### Note

**If you re-install IRsolution or install new version of it, it is strongly recommended that you uninstall the old software by the following procedure before installing the new software.**

1. Select the Control Panel on the Start menu of the Windows.
2. Select [Add / Remove Programs].
3. There are [IRsolution], [IRsolution x.xx] and [IRsolution Supplemental] on the window.
4. First, select [IRsolution Supplemental] then click the <Modify / Remove> button. Select [Remove] on the window then follow the prompts on the screen.
5. Next, select [IRsolution x.xx] then click the <Modify / Remove> button. Select [Automatic] on the window then follow the prompts on the screen.
6. Finally, select [IRsolution] then click the <Modify / Remove> button. Select [Remove] on the window then follow the prompts on the screen.
7. After uninstallation, close the [Add / Remove Programs] window and the Control Panel.