


### 7.7.3 [Hydrogen Leak Check] Screen

Press  (HOME) - [Peripherals] - [Hydrogen Leak Check Function] to display [Hydrogen Leak Check] screen.

Set this item when an optional hydrogen sensor (P/N: S221-88999-41) is connected.

▶▶ Reference "8.7.2 Hydrogen Sensor" P.319

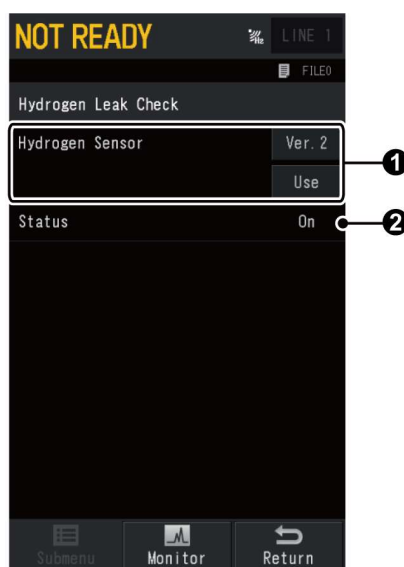

| No. | Item            | Description                                |                                                                                                                                                                            |
|-----|-----------------|--------------------------------------------|----------------------------------------------------------------------------------------------------------------------------------------------------------------------------|
| ①   | Hydrogen Sensor | Installation                               | Set [Ver.2] to display  (hydrogen sensor) icon on the screen, monitors hydrogen leak. |
|     |                 |                                            | Items N/A, Ver.1*1, Ver.2                                                                                                                                                  |
|     |                 |                                            | Default N/A                                                                                                                                                                |
|     |                 | Aspiration                                 | When you set [Use], aspirate atmosphere in the column oven.                                                                                                                |
|     |                 |                                            | Items Not Use, Use                                                                                                                                                         |
|     |                 |                                            | Default Use                                                                                                                                                                |
| ②   | Status*2        | Display the status of the hydrogen sensor. |                                                                                                                                                                            |

\*1 The [Ver.1] is not supported in the GC-2050.

\*2 The status is displayed when ① the hydrogen sensor is set to [Ver.2].