


# AOC-30 Series

If [AOC-30] is selected on the [Line Configuration] screen, the following screen appears.


[Line Configuration] Screen

■ **[Auto Injector] Screen**

To display [Auto Injector AOC-M], press  (HOME) - [AOC] after you install AOC-30 series on the instrument and configure AOC in the analytical line.

On this screen, you can check the status of AOC-30 series and set parameters for them.

Install AOC-30 series on the instrument and then configure AOC-M in the analytical line. For dual configuration, configure AOC-S (sub) in the analytical line, and switch lines to set up them.

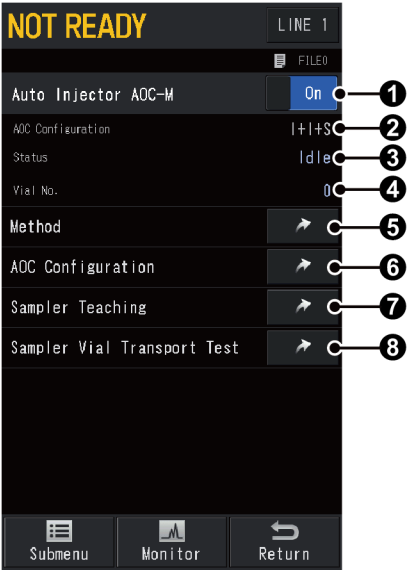

No.	Item	Description	
❶	Auto Injector Control	Select [On] to control AOC-30 series with the instrument.	
		Items	Off, On
		Default	Off