

[SCD Parts Counter] Screen

Press  (HOME) - [Detector] - [Submenu] - [SCD Parts Counter] to display [SCD Parts Counter] screen.
You can check usage time of SCD consumables.



No.	Item		Description	
①	Inner Pyrotube		Sets usage time of the inner pyrotube.	
		Usage Counter	Select [Yes] to display a message in the message display when usage time of the inner pyrotube exceeds the threshold.	
			Items	No, Yes
			Default	Yes
	Usage Time		Displays usage time of the inner pyrotube. Press [Reset] to reset usage time. When [Usage Counter] is set at [Yes] and usage time exceeds the threshold, the displayed color turns red.	
	Threshold	Sets usage time when you should replace the inner pyrotube.		
		Range	0 to 20000	
		Default	1000	
②	Outer Pyrotube		Sets usage time of the outer pyrotube.	
		Usage Counter	Select [Yes] to display a message in the message display when usage time of the outer pyrotube exceeds the threshold.	
			Items	No, Yes
			Default	Yes
	Usage Time		Displays usage time of the outer pyrotube. Press [Reset] to reset usage time. When [Usage Counter] is set at [Yes] and usage time exceeds the threshold, the displayed color turns red.	
	Threshold	Sets usage time when you should replace the outer pyrotube.		
		Range	0 to 20000	
		Default	1000	
③	SCD Ozone Scrubber		Sets usage time of the SCD ozone scrubber.	

	Usage Counter	Select [Yes] to display a message in the message display when usage time of the SCD ozone scrubber exceeds the threshold.			
		Items	No, Yes		
		Default	Yes		
	Usage Time	Displays usage time of the SCD ozone scrubber. Press [Reset] to reset usage time. When [Usage Counter] is set at [Yes] and usage time exceeds the threshold, the displayed color turns red.			
	Threshold	Sets usage time when you should replace the SCD ozone scrubber.			
		Range	0 to 20000		
		Default	8760		
	4	Vacuum Pump		Sets usage time of the vacuum pump.	
		Usage Counter	Select [Yes] to display a message in the message display when usage time of the vacuum pump exceeds the threshold.		
Items			No, Yes		
Default			Yes		
Usage Time		Displays usage time of the vacuum pump. Press [Reset] to reset usage time. When [Usage Counter] is set at [Yes] and usage time exceeds the threshold, the displayed color turns red.			
Threshold		Sets usage time when you should replace the vacuum pump.			
		Range	0 to 20000		
		Default	8760		