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A license under U.S. Patents 4,683,202, 4,683,195, 4,965,188, and 5,075,216 or their foreign counterparts, owned by Roche Molecular Systems, Inc. and F. Hoffmann-La Roche, Ltd. ("Roche"), has an up-front fee component and a running-royalty component. The purchase price of this product includes limited, nontransferable rights under the running-royalty component to use only this amount of the product to practice the polymerase chain reaction (PCR) and related processes described in said patents solely for the research and development activities of the purchaser when this product is used in conjunction with a thermal cycler whose use is covered by the up-front fee component. Rights to the up-front fee component must be obtained by the end user in order to have a complete license to use this product in the PCR process. These rights under the up-front fee component may be purchased from Applied Biosystems or obtained by purchasing an Authorized Thermal Cycler. No right to perform or offer commercial services of any kind using PCR including, without limitation, reporting the results of the purchaser's activities for a fee or other commercial consideration, is hereby granted by implication or estoppel. Further information on purchasing licenses to practice the PCR process may be obtained by contacting the Director of Licensing at Applied Biosystems, 850 Lincoln Centre Drive, Foster City, California 94404 or the Licensing Department, Roche Molecular Systems, Inc., 1145 Atlantic Avenue, Alameda, California 94501.

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