Supplemental (Ver. 1.00 SU2) Disk for AIMView32 Ver.1.00

Supplemental Disk (Ver1.00 SU2) for AIMView32 Ver1.00 (206-72418-91) is a supplemental program to make the AIMView32 software work on Windows 2000 Professional/XP Professional. Even the Restricted user accounts of Windows 2000 Professional/XP Professional can use AIMView32 by installing this program. Please install it when you control AIM-8800 with the computer of the Windows 2000 Professional/XP Professional environment. Please NEVER install this Supplemental disk when you use AIMView32 on Windows 95/98/98SE/Me environment.

Preparation for installation

- Confirm AIMView32 Ver.1.00 has been installed before it is executed.
- End AIMView32 when it is activated.
- NEVER install this Supplemental disk when you use AIMView32 on Windows 95/98/98SE/Me environment.

Installation

1. Please install the AIMView32 Ver. 1.00 software referring to a chapter installed of the manual of AIM-8800.
2. Insert the disk# 1 of this Supplemental Disk in the floppy disk drive.
3. Click the [Start] button of Windows, and then select the [Run] command.
4. Input following program onto the Edit box, and click the OK button.
A:¥Setup.exe
5. The setup wizard starts. Select the language of the installed program.
6. Follow the prompts of the wizard's instruction.

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