IMAGEREVEAL MS Installation Manual

1 Installing and Uninstalling

1.1 Installing

This section explains the procedure for installing all IMAGEREVEAL programs to your computer.

NOTE	 Note that once your IMAGEREVEAL is upgraded from Version 1.1 or earlier version to Version 1.2, it cannot be downgraded to Version 1.1 or earlier.
	 Depending on your anti-virus software, the installation may fail. In this case, disable your anti-virus software when installing this software.
	• When upgrading IMAGEREVEAL, be sure to exit the software before installing.
	 MATLAB Runtime compatible with your version of IMAGEREVEAL must be installed to ensure correct operation. After installing IMAGEREVEAL, do not uninstall only MATLAB Runtime.
	 When installing LabSolutions on the same computer, you must install IMAGEREVEAL before installing LabSolutions. The compatible combination of versions that can be installed on the same computer is as follows: IMAGEREVEAL Ver.1.20 or later LabSolutions LCMS Ver.5.109 or later LabSolutions Insight Ver.3.8 SP1 or later

Uninstall IMAGEREVEAL if the version is 1.1 or earlier.

Reference "1.2 Uninstalling IMAGEREVEAL"

NOTE Installing the latest version of IMAGEREVEAL while there is an existing version of IMAGEREVEAL (Version 1.1 or earlier) prevents you from launching IMAGEREVEAL. If Version 1.1 or an earlier version of IMAGEREVEAL is installed, uninstall it first before installing the latest version of IMAGEREVEAL.

Third-party trademarks and trade names may be used in this publication to refer to either the entities or their products/services, whether or not they are used with trademark symbol "TM" or "®".

© 2018-2022 Shimadzu Corporation. All rights reserved. Original version is approved in English.

Install MATLAB Runtime.

2

- 1 Copy the [MATLABRuntime] folder from the installation disk to any location on the PC where you want to install it.
 - NOTE MATLAB must be installed before installing IMAGEREVEAL. If the correct version of MATLAB Runtime is already installed, this procedure is unnecessary.
 If you unzip the [MATLABRuntime] folder without copying it to the PC where you are installing it, the unzip process will not complete successfully. Be sure to copy it. There is no need to copy it when installing with the iso file.
- 2 Select the copied [MATLABRuntime] folder and select [Extract All...] from the right-click menu.

The destination folder selection screen appears.

3 Click [Extract]. The extraction process is performed.

NOTE The extraction process may take several minutes to complete.

4 When the extraction is complete, double-click "setup.exe" in the extraction destination folder. Start the MATLAB installer.

NOTE It may take a few minutes for the installation window to appear.

5 Click [Next].

MATLAB Installation Window

6 Check the license agreement. If you agree, check the [Yes] checkbox and click [Next].

MATLAB License Agreement

7 Leave the installation folder as-is and click [Next].

MATLAB Folder Selection

NOTE Do not change the installation folder from the default.

8 Confirm the settings and click [Install].

MATLAB Confirmation Window

The installation starts.

📣 0% Complete			-		×
Installing					
	0%				
				Рац	JSE
		Cancel	∢ М	athWor	'ks•

MATLAB Installation Progress Window

9 After the installation ends successfully, click [Exit] to close the window.

MATLAB Installation Complete Window

3

Install IMAGEREVEAL.

1 Double-click "IMAGEREVEALSetup.exe".

NOTE If MATLAB Runtime is not installed, IMAGEREVEAL cannot be installed.

2 Click [Install].

Installation Window

The installation starts.

MAGEREVEAL Setup -	_		×
Setup Progress			
Processing: Microsoft .NET Framework 4.7.1			
	[Cance	el

Installation Progress Window

NOTE If the follo	owing window is displayed, click [Yes].	
	User Account Control × Do you want to allow this app to make changes to your device?	
	Setup Program Verified publisher: Shimadzu Corporation File origin: Hard drive on this computer Show more details	
	Yes No	
	[User Account Control] Window	

When the installation complete window is displayed, click [Restart] to restart your computer.

4

Installation Complete Window

NOTE • If you have installed anti-virus software that performs real-time scans when files are accessed, it may take more time to perform operations such as saving the project file or creating a data matrix. Configure the anti-virus settings so that the folders and files related to this software are exempt from the scans. • After this software is installed for the first time, you can use it as a 30-day demo version. However, the demo version has the following functional restrictions: "Demo" is displayed on MS images. • MS images, graphs, and data matrix tables cannot be exported. To use the software without these functional restrictions, or to continue using the software after 30 days, license authentication is required. When installing this software, the license management tool will be installed at the same time. Reference "2 License Authentication" P.9 • When using this software for the first time after installation, you must set the location to save the data matrix temporary file. Otherwise, a data matrix cannot be created.

1.2 Uninstalling IMAGEREVEAL NOTE Before uninstalling this software, make sure to exit the software. 1 In the [Start] menu, click [Settings] - [Apps]. 2 Click [IMAGEREVEAL], then click [Uninstall]. The uninstallation window is displayed. 3 Click [Uninstall].

Uninstallation Window

The uninstallation starts.

Uninstallation Progress Window

2

After the uninstallation ends successfully, click [Close] to close the uninstallation window.

Uninstallation Complete Window

1.3 Uninstalling the Software License Management Tool

NOTE If LabSolutions is installed, do not uninstall the Software License Management Tool. Doing so prevents the license authentication for LabSolutions.

1 In the [Start] menu, click [Settings] - [Apps].

Click [ActivationModulePackage], then click [Uninstall].

NOTE If the following	window is displayed, click [Yes].	
	User Account Control Do you want to allow this app to make changes to your device?	×
	9a9f8.msi Verified publisher: Shimadzu Corporation File origin: Hard drive on this computer Show more details	
	Yes No	
	[User Account Control] Window	

The uninstallation window is displayed, and the uninstallation process is completed.

2 License Authentication

Authenticate your license by following the instructions on authenticating the software license for the Software License Management Tool.

- NOTE If you have a time-limited license, note that it cannot be canceled during its term. Irrespective of whether you actually use the software, once your license is activated, it will expire after its valid term starting from the date of activation.
 - When installing IMAGEREVEAL and LabSolutions on the same computer, you will see the message "Undefined license included." in the LabSolutions' License Information window. This does not create any problem for correct application operation.

Follow the steps below to display the software license authentication procedure manual.

- From the [Start] menu, select [All Apps List] [Shimadzu License Management Tools] - [Software License Management Tool].
 If the [User account management] window is displayed, click [Yes].
- 2 Click the [Help] button in the Software License Management Tool window. The software license authentication procedure manual appears.

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▲ MATLAB Runtime インストーラー	-		×
このアプリケーションは MATLAB Runtime をインストールします。			
	MA	TLA RUNTIME R2019a	AB⁺
			L.
MATLAB and Simulink are registered trademarks of The MathWorks, Inc. Please see mathworks.com/trademarks for a list of additional trademarks. Other product or brand names may be trademarks or registered trademarks of their respective holders.			
WARNING: This program is protected by copyright law and international treaties. Copyright 1984-2019, The MathWorks, Inc. Protected by U.S. and other patents. See MathWorks.com/patents			
< 戻る 次へ > キャンセル	📣 M	athWor	ks•

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5 单击 [Next]。

A MATLAB Runtime Installer	-		×
This application will install MATLAB Runtime.			
	MA	TLA RUNTIME R2019a	AB*
			L.
MATLAB and Simulink are registered trademarks of The MathWorks, Inc. Please see mathworks.com/trademarks for a list of additional trademarks. Other product or brand names may be trademarks or registered trademarks or their respective holders.			
WARNING: This program is protected by copyright law and international treaties. Copyright 1984-2019, The MathWorks, Inc. Protected by U.S. and other patents. See MathWorks.com/patents			
< <u>R</u> ack Next > Cancel	📣 М	athWor	'ks•

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📣 0% Complete			-		×
Installing					
	0%				
				Pau	ise
		Cancel	- 4 N	(athMar	101

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📣 Installation Complete	-		×
Installation is complete.			
	M	RUNTIME R2019a	AB*
			L.
< Back Figish Cancel	- 4 M	fathWor	'ks•

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✓ 过释 如果显示以	下窗口,请单击[Yes]。	
	User Account Control Do you want to allow this app to make changes to your device?	×
	Verified publisher: Shimadzu Corporation File origin: Hard drive on this computer Show more details Yes	
	用户帐户控制窗口	

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2 单击 [ActivationModulePackage] 后,单击 [卸载]。

♥ 注释	显示以下窗口时,	请单击[是]。	
		User Account Control X Do you want to allow this app to make changes to your device?	<
		9a9f8.msi Verified publisher: Shimadzu Corporation File origin: Hard drive on this computer Show more details Yes No	
		用户帐户控制窗口	

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本文中に記載されている会社名、製品名、サービスマーク、およびロゴは、各社の商標および登録商標 です。本文中では「TM」、「®」を明記していない場合があります。 © 2018-2022 Shimadzu Corporation. All rights reserved.
在本说明书中记载的公司名称,产品名称,服务标志及logo是各公司的商标及注册商标。另外,本 说明书中不对[™]、®标记做明确说明。 本说明书是日文版《IMAGEREVEAL MS安装手册》(225-28894)(修订版C 2021年7月)的译文。 © 2018-2022 Shimadzu Corporation. All rights reserved.