

BX-K、BW-K シリーズをこれからお使いの皆様へ

このたびは BX-K、BW-K シリーズをお買い上げいただき誠にありがとうございます。

最近、応答性を重視した設定をご希望されるお客様が多くいらっしゃいます。それにともない、当社工場出荷時に【平均化処理の設定】(取説 P.26)を本来の「自動モード 22」から、「はかりとりモード 23」(モード変更方法は取説 P.14~16)にあらかじめ設定変更してあります。

もし設置環境の外乱が大きく測り難い場合や、はかりとり以外の作業にて本製品をご使用の場合は取扱説明書の 2.8.2 項【平均化処理の設定】(取説 P.26)をご参照いただき、本製品の表示の安定性と応答性の程度を適切に設定してください。なおメニューリセット(取説 P.19)を行うと「自動モード 22」に平均化処理の設定が自動的に行なわれますので、ご注意ください。

Dear customers,

Thank you for purchasing BX-K/BW-K series.

In order to have good weighing data quickly, we changed the factory preset of “Averaging” to “POURing” (Menu Map item [22]) from “AUTOMatic” (Menu map item [23]) in Environment Menu. (Refer “2.8.2 Averaging”, P.30 of the instruction manual.)

* Also refer P13-20 for menu map.

If the balances is installed in an unstable environment or if you need higher stability, please change the setting appropriately depending on the environment or the usage.

Note: If you use the menu reset function (“2.4.5 Returning to the Default Settings”, P.22), it will be set as “AUTOMatic”