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RDB-200 Series Remote Display Unit Dedicated for Shimadzu Electronic Balances Instruction Manual

1. Introduction

The RDB-200 series remote display unit is connected to a Shimadzu electronic balance and is used to read the displayed contents and operate the balance from a location away from the balance unit.

! Caution Do not connect the display unit to equipment other than compatible Shimadzu electronic balances.
Connecting it to other equipment may cause a failure.

2. Installation

Do not use the product in places such as those listed below.

- Places with dust, vibration, electromagnetic waves, or magnetic fields
- Places exposed to direct sunlight and places near heat sources such as a heater.
- Places with a drifting corrosive gas
- Places with an extremely high or low temperature or high or low humidity

3. Connecting with a Balance and Removing

Connecting procedure

- (1) Pull out the AC adapter of the balance main unit from the power supply outlet.
- (2) Properly connect the connector of the RDB-200 series unit to the balance main unit.
- (3) Insert the AC adapter into the power supply outlet and turn on the power.
- (4) Press the [POWER] key on the balance.

Removing procedure

- (1) Press the [POWER] key on the balance to put the balance into standby.
- (2) Pull out the AC adapter of the balance main unit from the power supply outlet.
- (3) Remove the connector of the RDB-200 series unit from the balance main unit.

! Caution Do not insert or remove the connector of the RDB-200 series unit with the power of the balance main unit turned on.
Incorrect operation may cause a failure.

4. Key Operations (RDB-201 only)

Key	Operation Function
[POWER]:	Turn on/off the power, bypass menus
[MENU]:	Select menus, calibrate the balance
[TARE]:	Tare, determine menus
[RANGE]:	Change the measurement range (This key functions as the [TARE] key for models other than applicable models.)
[UNIT]:	Change the unit
[PRINT]:	Print data using a printer

Key operation functions of the RDB-201 are the same as those of the balance main unit to which the unit is connected. For detailed functions of each key, refer to the instruction manual provided with the balance main unit.

5. Specifications

RDB-201	Remote display unit, with operation keys
RDB-202	Remote display unit only
Length of connection cable	Standard approx. 2 m
Housing size	Approx. W165 × D100 × H50 mm
Weight	Approx. 600 g

6. Peripheral Equipment

(1) Angle adjuster and wall hook set (P/N: 321-55953)

Mounting this product on the RDB main unit enables you to adjust the angle of the display unit. It also enables you to hang the display unit on the wall.

For detailed installation procedure and how to use it, refer to the instruction manual provided with the angle adjuster and wall hook set.

(2) Stand for remote display (P/N: 321-55954-01) (1 m high)

This stand is dedicated for the RDB-200 series units.

It improves the visibility and operability of the display unit in a workspace with no desk or table. The angle of the display unit can be adjusted.

For detailed installation procedure and how to use it, refer to the instruction manual provided with the stand for remote display.