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# High Performance Packed Column for High Performance Liquid Chromatograph

Shim-pack

# SCR-102H

## Instruction Manual

### ■ Introduction

The Shimadzu Shim-pack SCR-102H is a column of 8 mm *i.d.* 30 cm long packed with H type cation exchanger consisting of semi-rigid styrene-divinylbenzene copolymer.

Shim-pack SCR-102H is a especially suitable column for the analysis of the organic acids based on ion exclusion chromatography.

### ■ Column Performance


Column test report and chromatogram are included. Guaranteed efficiency in terms of theoretical plate numbers is 15,000.

### ■ Column Installation

- For installing the column with stainless steel piping, the following connection parts are required.

Item Name	P/N	Comments
Male nut 1.6 MN	228-16001	2/pkg
Ferrule 1.6 F	228-16000	2/pkg

- Connect as in the Figure.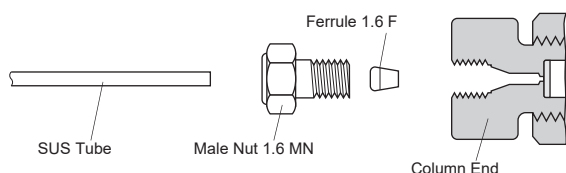

Do not tighten the joints excessively !

### ■ Mobile Phase Solvent

- As a mobile phase solvent, 0.1 % of phosphoric acid or perchloric acid aq. solution can be used.
- Some samples which tend to have hydrophobic property retain longer due to adsorptive interaction. The addition of Ethanol to mobile phase will bring faster retention. (The addition of Ethanol should be less than 20 %.)

### ■ Column Handling Precautions

- Mobile phase flow in the direction indicated on the column label only.
- Maximum usable pressure is 50 kg/cm<sup>2</sup>. Higher pressure will degrade the column performance. Do not allow high pressure by delivering the solvents very fast.

- Appropriate flow rate is 1.5 mL/min under the temperature of 40-85 °C. When used at lower than 40 °C, pay attention the flow rate to keep the maximum usable pressure. (Particularly, in cold season.)
- When increase or decrease the column temperature, set the flow rate at 0.2 mL/min until the column temperature reaches at settled value.
- Avoid the vigorous change of pressure and flow rate to protect the column bed.
- Keep the maximum usable column temperature of 85 °C.
- Dust particles in mobile phase of sample solution lead to plugging at the column inlet filter or at the top of column bed. Filter the mobile phase and sample solution prior to use, if necessary.
- Metal ions and amino compounds etc. in sample or mobile phase solution may degrade the separability of columns in which cation exchange resin is packed. For guard from such trouble, use exclusive column [Guard Column SCR-102H, part number 228-17924-91].
- The solvent filled in the column at the time of shipment is 0.1% Phosphoric acid aqueous solution.
- Replace the enclosed mobile phase within 3 months.
- Before using a new or unused column for more than 3 months, flush the column overnight with the mobile phase by setting the flow rate to approximately 0.1 mL/min.

### ■ Column Storage

- For long term storage wash the column and fill it with mobile phase solvent followed by plugs to seal the column.

### ■ Technical Support

It is the customer's responsibility to develop and validate analytical conditions for a particular application. However, Shimadzu offers technical support by e-mail and phone for customers who need help.

Write specific questions to [analytic@group.shimadzu.co.jp](mailto:analytic@group.shimadzu.co.jp) or call your local representative.

\*The contents of this instruction sheet are subject to change without notice.