Shimadzu Packed Column for HPLC

Shim-pack™ SCR-101C

Instruction Manual

■ Introduction

The Shimadzu Shim-pack SCR-101C is a column of 7.9 mm I.D. and 300 mm long stainless steel tube packed with Ca type cation exchanger consist of styrenedivinylbenzene copolymer.

Shim-pack SCR-101C is an exclusive use column for the analysis of sugars, glycitols and polyalcohols based on gel chromatography or ligand exchange chromatography.

■ Operating Care

- Mobile phase flow in the direction indicated on the column label only.
- Maximum usable pressure is 4 MPa. Higher pressure will degrade the column performance. Do not allow high pressure by delivering the solvents very fast.
- Appropriate flow rate is less than 1 mL/min under the temperature of 40 to 85 °C
- Avoid the vigorous change of pressure and flow rate to protect the column bed.
- Column temperature of 80 °C brings better separation. Keep the maximum usable temperature of 90 °C.
- When increasing or decreasing the column temperature, set the flow rate at 0.2 mL/min until the column temperature reaches at set value.
- Do not subject to shock such as from dropping, etc.
- Dust particles in mobile phase or sample solution lead to plugging at the column inlet filter or at the top of column bed. Filter the mobile phase and sample solution prior to use, if necessary.
- Metal ions and amino compounds etc. in sample or mobile phase solution may degrade the separation of columns in which cation exchange resin is packed.
For guard from such trouble, use exclusive column [Guard Column SCR(C), part number 228-17891-91].
- The solvent filled in the column at the time of shipment is pure water.

■ Selection of Mobile Phase

- As a mobile phase solvent only pure water can be available.

■ Maintenance of Column


- For long term storage fill it with pure water followed by plugs to seal the column.
- If the separation became worth, wash the column with 0.1 M Calcium Nitrate at the flow rate of 0.5 mL/min, 80 °C for 3 hours. Then wash with pure water for 1 hours. If the column did not recover, then considered the life is end.

■ Column Connection

- Tubing kit available for column connection is listed below.

Item Name	P/N	Comments
Ferrule 1.6F 316L	228-16000-10	1/pkg
Male nut 1.6MN	228-16001	1/pkg

- Observe the appropriate connection method as illustrated in the Figure.

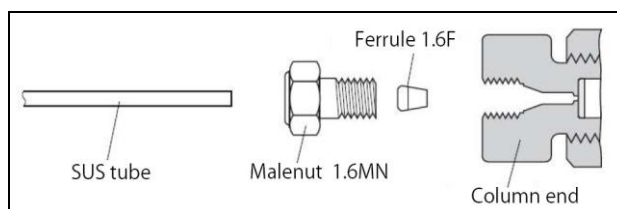- Do not overtighten as this will result in damage to the fittings.

■ Technical support

Shim-pack SCR columns are manufactured, inspected, packaged and shipped under strict standards of quality control. Should you find any defect in performance, please contact your local Shimadzu representative, who will ensure your complete satisfaction.

We regret that we cannot guarantee the lifetime of columns, also that we cannot accept any claim when performance has deteriorated due to noncompliance with the operation procedures elucidated above, or as a result of normal aging.

※ The contents of this instruction sheet are subject to change without notice.

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