## Shimadzu Packed Column for HPLC

# Shim-pack™ SCR-101P

## Instruction Manual

### ■ Introduction

The Shimadzu Shim-pack SCR-101P is a column of 7.9 mm I.D. and 300 mm length stainless steel tube packed with Pb type cation exchanger consisting of styrene-divinylbenzene copolymer.

Shim-pack SCR-101P is suitable for the analysis of sugars, sugar alcohols and polyalcohols by gel chromatography and ligand exchange chromatography.

### ■ Column handling precautions

- The flow direction is shown with an arrow on the column label (→). When installing the column, ensure that the mobile phase flow direction matches the direction on the column.
- Maximum usable pressure is 4 MPa. Extreme pressure may cause the degradation of the column performance. Do not exceed an operating pressure of 4 MPa.
- The recommended flow rate is less than 1 mL/min between 40 to 85 °C.
- Avoid excessive pressure fluctuation. This increases pressure rapidly at the column inlet, which may cause premature column deterioration.
- The recommended operating temperature of the Column is 80 °C for separating the compounds. The maximum operating temperature is 90 °C.
- When increasing or decreasing the column temperature, set the flow rate at 0.2 mL/min until the column temperature reaches at the set value.
- Do not drop or bump the column to avoid damage to the column.
- The most common cause of increased column backpressure or split peaks is clogging of the inlet frit by sample particulates. Filter the mobile phase and sample solution before flushing the flow line and connecting the column, if necessary.
- The compounds such as metal ions and amino compounds in sample solution and mobile phase may degrade the separation of columns because cation exchange resin is packed into the column. Installing a guard column can prevent also prevent column clogging and prolong lifetime. The part number of the Guard Column SCR(P) is 228-17892-91].
- The shipping solvent is 100% water.

### ■ Mobile phase selection

- Water can be used as the usable mobile phase.
- Do not add any organic solvents to the mobile phase.

### ■ Cleaning and storing the column


- When storing the column, fill the column with water, securely plug both ends and store the column at ambient temperature.
- When the low-resolution chromatography is observed wash the column with 0.1 M Lead Nitrate solution at the flow rate of 0.5 mL/min, 80 °C for 3 hours. Then wash the column with water for 1 hours. If the flushing procedure does not solve the problem, replace the column.

### ■ Column Connection

- The following parts are required for column connection.

| Item Name         | P/N          | Comments |
|-------------------|--------------|----------|
| Ferrule 1.6F 316L | 228-16000-10 | 1/pkg    |
| Male nut 1.6MN    | 228-16001    | 1/pkg    |

- Connect the column as shown in the figure below.

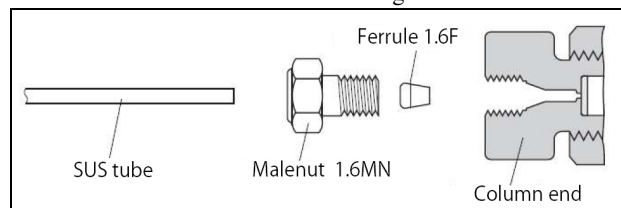- Do not overtighten the nut. It may cause damage to the column inlet fitting.

### ■ Technical support

Shim-pack SCR series columns are manufactured, inspected, packaged and shipped under strict standards of quality control. Shimadzu offers technical support for customers who need to help. Contact your local representative or a nearby Shimadzu sales office if needed.

※ The contents of this instruction sheet are subject to change without notice.

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