High Performance Packed Column for High Performance Liquid Chromatograph

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SCR-102H

Instruction Manual

- Shim-pack Guard column SCR-102H is a guard column (6 mm i.d., 5 cm long) packed with H ion substituted exchanger consist of sulphonated styrene-divinylbenzene copolymer.
- Guard column SCR-102H is exclusively used in combination with Shim-pack SCR-102H separation column. Refer the manual of Shim-pack SCR-102H column as well.
- Tubing kit available for column connection is listed in Table 1.

Table 1

Item Name	P/N	Comment
Column connector kit ASSY	228-17943-92	Ferrule 1.6 F × 2 Male nut 1.6 MN × 2

- Observe the appropriate connection method as illustrated in Fig. 1.

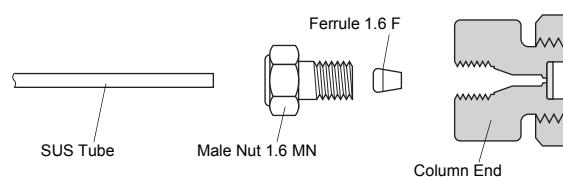
Fig.1 : Column Connection

Do not overtighten as this will result in damage to the fittings.

■ Technical Support

It is the customer's responsibility to develop and validate analytical conditions for a particular application. However, Shimadzu offers technical support by e-mail and phone for customers who need help.

Write specific questions to analytic@group.shimadzu.co.jp or call your local representative.

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*The contents of this instruction sheet are subject to change without notice.

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 **SHIMADZU CORPORATION**

ANALYTICAL & MEASURING INSTRUMENTS DIVISION

1, Nishinokyo-Kuwabara-cho, Nakagyo-ku, Kyoto, 604-8511, Japan