High Performance Packed Column

for High Performance Liquid Chromatograph

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
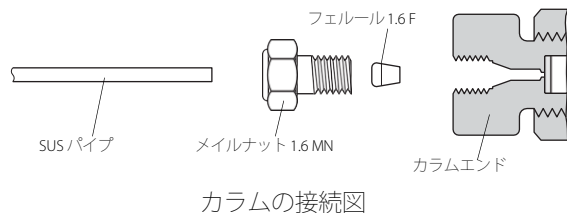
Shim-pack Guard column SCR(C)

Instruction Manual

- Shim-pack Guard column SCR(C) is a guard column(4 mm I.D., 50 mm L.) packed with Ca ion substituted cation exchanger consisting of styrene-diviny-benzene copolymer.
- Guard column SCR(C) is exclusively used in combination with Shim-pack SCR-101C. Refer to the manual of Shim-pack SCR-101C column as well.
- The tubing kit available for column connection is listed below.

Item Name	P/N	Comment
Column connector kit	228-17943-92	Ferrule 1.6 F x 2 Male nut 1.6 MN x 2

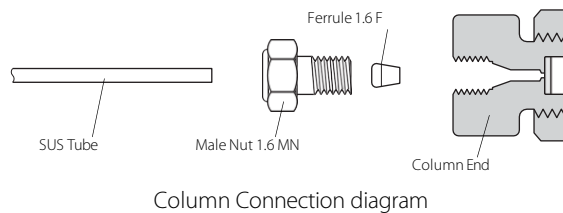- When installing the guard column, please refer to the diagram below to connect the joints to both ends of the column. Be careful not to over-tighten, as excessive tightening may lead to damage at the connection part.

■ Disposal Precautions

Separate the column from general industrial or household waste and dispose of it in accordance with local government regulations.

■ Technical Support

If you encounter any performance defects, please contact your local Shimadzu sales office for assistance. However, please note that we cannot guarantee the lifetime of the columns, nor can we accept any claims when performance has degraded due to non-compliance with the operation procedures outlined above, or as a result of normal aging.

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*Please note that the appearance, specifications, and the contents of this instruction sheet are subject to change without notice.