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## EP-100(321-78181) EP-110(321-78182)

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## Corrections to Instruction Manual for EP-100(321-78183), EP-110(321-78184)

The following is an addition to "6. Trouble shooting" in the EP-100 instruction manual.

Symptom	Cause	Remedy
The date and time have changed.	The built-in clock backup battery is low.	Plug in the AC adapter and turn it on more than 8 hours.
	Unintended settings were made. Press  and  simultaneously to print the setting change history. The cause of the change will be identified and printed as follows:  <b>User:</b> The date and time were changed by user operating. <b>CMD_USB:</b> The date and time were changed by EasySetup program. <b>POWER LOST:</b> The date and time were by the RTC backup battery power failure.	Set the correct date and time.
RTC ERROR	Internal clock data error	Turn the power switch back on.

The following is an addition to "7. Trouble shooting" in the EP-110 instruction manual.

Symptom	Cause	Remedy
The date and time have changed.	The built-in clock backup battery is low.	Plug in the AC adapter and turn it on more than 8 hours.
	Unintended settings were made. Press  and  simultaneously to check the setting change history. Press  to print the change history. The cause of the change will be identified and displayed as follows:  <b>User:</b> The date and time were changed by user operating. <b>CMD_USB:</b> The date and time were changed by EasySetup program. <b>POWER LOST:</b> The date and time were by the RTC backup battery power failure.	Set the correct date and time.
RTC ERROR	Internal clock data error	Turn the power switch back on.

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