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The screenshot displays the IMAGEREVEAL software interface. The main window is titled "差異解析" (Difference Analysis) and shows a "MSイメージ一覧" (MS Image List) panel on the right. This panel contains a grid of MS images, with the first image, "Testicle_9AA_PL_SL_5x_1_AREA01.mdx", highlighted by a yellow box and labeled "選択" (Select). The interface also features a "グラフ" (Graph) section showing a mass spectrum plot with peaks labeled at m/z 795.485, 795.521, 796.524, and 797.524. The "ROIリスト" (ROI List) table is visible at the top, listing various ROIs with their m/z values and PCA coordinates.

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The screenshot displays the IMAGEREVEAL software interface. The main window shows a data table with columns for No., 使用, タグ, ラベル, m/z, PCA-水平軸, PCA-垂直軸, Group A-平均, and Group B-平均. The selected m/z is 795.48492. Below the table, a box plot shows the intensity distribution for Group A (red), Group B (blue), and Group C (green). The y-axis is labeled 'Intensity' and ranges from 0 to 1,600,000. Below the box plot, two histograms are shown: one for Group A (red bars) and one for Group B (blue bars). The y-axis for the histograms is labeled 'Count'.

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